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FUNCTION

TOF emergency showers and eye washers comply with CE - DIN 12899 - UNI 9608 standards and they feature a high water flow, but a soft water spout for an effective decontamination of injured body parts, which have accidentally been hit by acid or corrosive chemicals or another harmful substance. Eye washers are indispensable in all areas at risk. They are easy to recognize by their green or red colour and a sign with a pictogram sign according to the standards.

INSTALLATION AND OPERATING INSTRUCTIONS

Floor installation in all areas at risk. The valve of the eye washer can be opened either by pushing down the pedal or the square hand lever with the identification sign. The shower head valve is opened by pulling the triangular tie rod.

The valves stay open during the whole flushing process.

TECHNICAL DETAILS OF CONNECTION

G1/2 - G1 (according to the type) to be connected to the drinking water-main.

OUTLET: G1

MINIMUM WATER PRESSURE: 2 bar

WATER FLOW

12 l/min for eye-washers (according to the standards).

30 l/min for showers (according to the standards).

WATER TEMPERATURE: 15°C - 35°C

Warning: excessive water temperatures might accelerate some chemical reactions. In this case we suggest to ask a specialist for advice.

MAINTENANCE

According to local standards of the country where the showers and eye washers are located.

REPAIRS

Emergency showers and eye washers may only be repaired in our premises or by skilled and qualified staff. TOF does not undertake any responsibility for alterations or repairs without its consent.

WARRANTY

TOF gives warranty in accordance with the sales contract and the general sales conditions. All products carry a warranty on assembly and manufacturing faults during six months after the delivery date indicated on the shipping documents. The warranty is strictly limited to faults discovered under normal circumstances and correct use. No warranty claim will be accepted for defects caused by alterations or incorrect use by the customer.

IMPORTANT

These instructions can be applied to all TOF showers and eye washers of series 1100. The drawings do not always show the supplied item.

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